

# SMART system<sup>5</sup>

File Name : Food\Fat\Mayo.doc

Revision 8/1999

## Application Note


Sample Type: **Mayonnaise**

Sample Prep Required: Stir the sample well.

Parameters: Sample Size: 2-3 grams

	<u>Moisture</u>	<u>Fat</u>
Power:	100 %	100 %
Delta Weight*:	0.5 mg	0.5 mg
Delta Time*:	10 seconds	10 seconds
Max Time:	3 minutes	2 minutes
Max Temp:	100 °C	

Spread Technique: Spread the sample quickly and smoothly over 90% of the pad surface area.

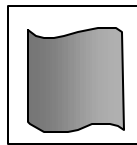

Notes: \*The above parameters are suggested starting points and may need to be optimized to meet certain product specifications.

✧ For General Guidelines and Operation Consult the SMART System<sup>5</sup> Operations Manual.

**CEM Corporation**  
3100 Smith Farm Road  
Matthews NC 28104