


# SF<sub>6</sub> Purity Analysis by Micro GC

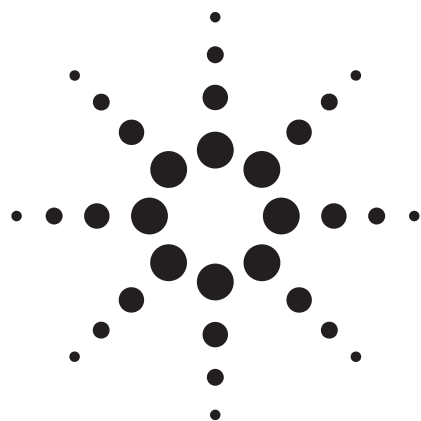

Application

Specialty Chemical

## Author

Wenmin Liu  
Agilent Technologies, Inc.  
412 Ying Lun Road  
Waigaoqiao Free Trade Zone  
Shanghai 200131  
China

## Abstract

**An SF<sub>6</sub> purity analysis method is developed in this application. PLOT Q and GasPro columns are suitable for the analysis of contaminants in SF<sub>6</sub>. On the GasPro channel, sulfur compounds are detected and the total analysis time is less than 2 minutes.**

## Introduction

SF<sub>6</sub> circuit breakers operate to switch electric circuits and equipment in and out of the system. These circuit breakers are filled with compressed sulfur-hexafluoride gas, which is a dielectric used in high-voltage electrical equipment as an insulator and/or arc-quenching medium; the gas also interrupts the current flow when the contacts are open. SF<sub>6</sub> gas is expensive and has been designated as a greenhouse gas by the EPA.

A reliable field method is important to assess whether reclaimed SF<sub>6</sub> is acceptable for continued use as in-service gas in high-voltage equipment, thus preventing failures, extending equipment life, reducing maintenance costs, and increasing personnel safety.

In this application, a method based on the Agilent 3000 Micro GC was developed to accurately determine concentrations of impurities and decomposition products in SF<sub>6</sub> to ensure reliable operation of electrical equipment.

## Experimental

An Agilent 3000 Micro GC with two-channel configuration: PLOT Q (8 m × 0.32 mm) and GasPro (10 m × 0.32 mm) was used.

The SF<sub>6</sub> sample was provided by Hebei Electric Power Research Institute.

The chromatographic conditions for the two channels are shown in Table 1.

**Table 1. Analytical Conditions of Micro GC**

Channel	PLOT Q	GasPro
Sample inlet temperature (°C)	80	80
Injector temperature (°C)	80	80
Column temperature (°C)	60	60
Inject time (ms)	100	100
Column pressure (psi)	20	20

## Results

The PLOT Q channel and the GasPro channel are both suitable for the purity analysis of SF<sub>6</sub>. Figures 1a and 1b show the chromatograms obtained from these two columns. Figure 1c is a zoomed-out window of Figure 1b from the time range of 1.0 to 2.6 minutes. An unknown sulfur




Agilent Technologies

compound is detected in the GasPro channel. The total analysis time is less than 2 minutes, providing excellent cycle time. The peak area repeatabilities (shown by RSD, n = 6) of these two channels are shown in Table 2. The detection limits of this method are approximately 20 ppm, which meets the user requirement for contaminant analysis in SF<sub>6</sub>.

**Table 2. The Repeatabilities of Two Channels**

RSD (n = 5)	Air (%)	CF <sub>4</sub> (%)	SF <sub>6</sub> (%)	CO <sub>2</sub> (%)
PLOT Q	2.80	0.19	0.16	0.98
GasPro	3.19	0.68	0.12	

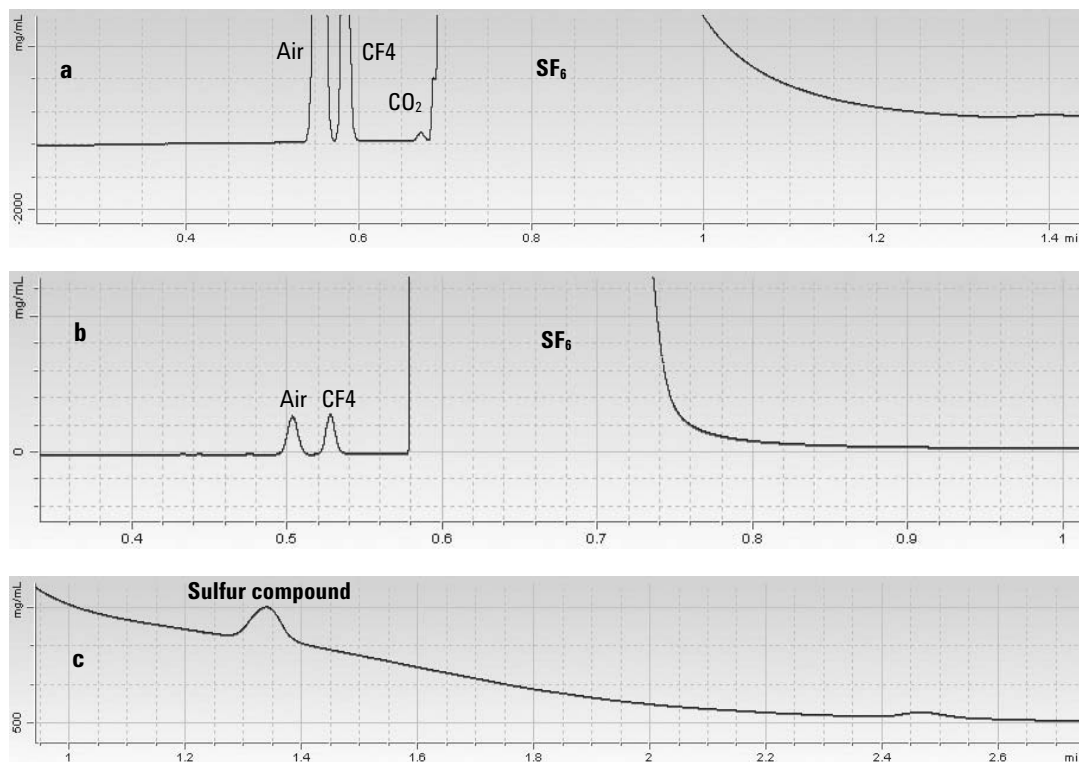

**Figure 1. Chromatograms in PLOT Q (a) and GasPro (b and c) channels.**

## For More Information

For more information on our products and services, visit our Web site at [www.agilent.com/chem](http://www.agilent.com/chem).

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc. 2007

Printed in the USA  
September 21, 2007  
5989-7430EN